




MPMS Application Note 1014-819


## Connecting a Straw Adapter to the Standard Transport Rod

1. Slide the straw adapter (Part #4084-234) onto the standard transport rod until the holes in the adapter are aligned with the holes in the end of the standard transport rod.

**Note:** The straw adapter's internal threads don't match the threads of the standard transport rod. Disregard the straw adapter's threads since the inner diameter of the straw adapter is large enough to accommodate the standard transport rod.

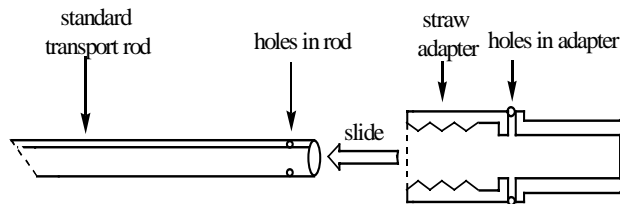

2. Connect the straw adapter to the standard transport rod by inserting a piece of copper buss wire through the aligned holes.

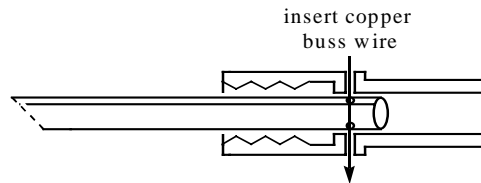

3. Once the straw adapter is secure to the rod, place a straw on the straw adapter.

